

# Data Sheet | Specifications

# **Tensormeter Model RTM1**

# Sheet and Hall resistance measurements: automated, ultra-precise, offset-free



Tensormeter model RTM1 front panel

The Tensormeter is designed for automated precision measurements of resistances and voltages. It unites the benefits of Lock-in Amplifiers and Source/Measure Units through an innovative flexible architecture based on an integrated matrix switch. Tensormeter RTM1 enables the automated recording of the complete Resistivity Tensor ( $R_x$ ,  $R_y$ ,  $R_H$ ) with one single device, even on unpattern thin films. With excellent AC and DC performance, it covers the range from Nano-Ohm to Giga-Ohms with at least 8 digits of dynamic range.

### **Application fields**

Materials research and characterization

- solid state physics
- semiconductor physics
- magnetism
- flexible electronics
- spintronics
- new functional electronic materials and devices

#### Industrial R&D and wafer/device testing

- microelectronic devices
- memory devices
- transistors, diodes
- LED/OLED
- solar cells
- displays, TCO
- sensors






#### **Benefits:**

- Replaces all standard devices for electrical characterization measurements (e.g. Lock-in Amplifier, SMU, DMM).
- Overcomes the limitations of stationary 4-point measurements by an integrated Matrix Switch.
- Offers presets for van-der-Pauw and Resistance Tensor measurements and allows for full user configurability.
- Makes complex sample preparation unnecessary (e.g. lithographic structuring).
- Allows for easy connectivity to many different measurement setups (e.g. probe stations, cryostats, vacuum systems).
- Saves measuring time and enhances sample throughput.

#### **Features:**

- Reconfigurable device architecture based on an integrated switching matrix
- 8 user-defined channels (BNC connectors), whose function (input or output) can be freely determined
- Conventional AC and DC 4-wire measurements with fixed connections (Kelvin/ Hall geometry)
- AC and DC measurements with alternating connections (van-der-Pauw geometry) with one device
- Simultaneous measurement of exactly separated absolute values for longitudinal and transverse resistances without lithographic patterning
- Software presets for common measurement modes, but any user-specific switching sequences can be specified
- TCP-based communication, easy integration in any environment (e.g. Labview, C, Python)

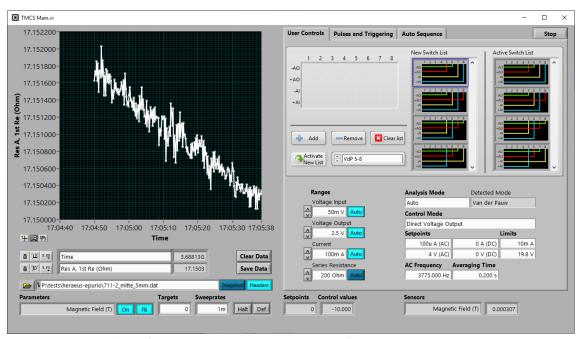

Graphical user interface of proprietary Tensormeter control software






## **Typical measurement examples:**

- Ultra-low noise and high stability AC & DC 4-wire measurements in standard geometries (Kelvin and Hall layouts)
- Van-der-Pauw switched connection 4-wire measurements on irregular, unstructured thinfilm samples
- Zero-Offset Hall 4-wire measurements (exact separation of longitudinal and transverse resistance even with unstructured samples)
- Sub-ppm relative resistance change measurements
- Ratiometric resistance measurements to eliminate sample and device drifts
- High drive harmonic distortion measurements, pulse & measure routines

#### Low Resistive Sensors and Specimen


Differential Input Noise Spectrum of a resistive sensor. Ultra-low wideband & 1/f noise AC measurements allow accurate sensor characterization and operation.

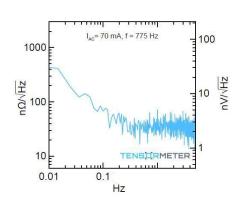


#### Zero-Offset Hall: Eliminate Drift and Parasitics

Differential Input Noise Spectrum of a Hall measurement on a thin film sample. The Zero-Offset Hall preset of the Tensormeter eliminates thermal drift and allows long integration and orders of magnitude improved sensitivity compared to regular 4-wire Hall measurements.

Loss of magnetization during warmup of an antiferromagnetic sample monitored in Hall Resistance. The Zero-Offset Hall preset of the RTM1 (top) clearly shows the loss of signal. On the contrary, parasitic signal contributions overshadow the useful magnetization signal in a regular 4-wire Hall measurement of the same sample (bottom).

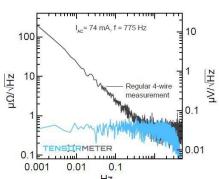

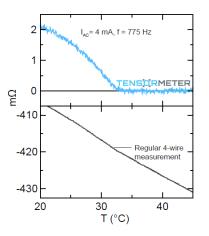





# **Electrical Specifications (typ.)**

#### **Outputs**

- Symmetrical Output: DC to 20 kHz, ±20 V, ±100 mA
- Output noise: < -140 dBFS
- Pulse / arbitrary function output with 10 μs resolution

#### Switch Matrix

- Fully controllable integrated 8x4 switching matrix
- Maximum switching frequency: 100 Hz
- LEDs indicate the switch configuration (can be dimmed or switched off)

#### Measurements

- Default FET option input voltage noise:  $5 \text{ nV}/\sqrt{\text{Hz}}$ ,  $1 \text{ fA}/\sqrt{\text{Hz}}$
- Optional BJT replacement input voltage noise:  $3 \text{ nV}/\sqrt{\text{Hz}}$ ,  $400 \text{ fA}/\sqrt{\text{Hz}}$
- Resistance measurement noise floor:  $50 \text{ n}\Omega/\sqrt{\text{Hz}}$  (direct BJT input)
- Resistance measurement noise floor:  $10 \text{ n}\Omega/\sqrt{\text{Hz}}$  (Transformer coupled FET input)
- Gain change with temperature: 100 ppm/K
- Gain change with temperature in ratiometric mode: <1 ppm/K
- DC offset voltage change with temperature: ±1 μV/K, ±3 ppm/K
- DC offset voltage change with temperature in differential mode: <0.1 ppm/K</li>

#### Ranges

- 4 User/Auto selectable Series Resistors from 200  $\Omega$  to 200  $k\Omega$
- 4 User /Auto selectable Output Voltage Measurement Ranges from 2.5 V to 20 V
- 4 User /Auto selectable Current Measurement Ranges from 100 nA to 100 mA
- 8 User /Auto selectable Input Voltage Measurement Ranges from 100 mV to 20 V
- Base Precision within Range: <1 ppm of full range for Output Voltage and Current</li>
- Base Precision within Range: <0.1 ppm of full range for Input Voltage
- Continuous dynamic range within one Range: > 8 digits

#### <u>Other</u>

- Arbitrary function reference input/output: single-ended ±10 V
- Trigger input/output: single-ended 5V TTL





# **Hardware/software specifications**

- 19" rack-mountable device, 3 height units, 25 cm depth
- TCP-based user connection
- Windows driver and client communication examples for LabView and Python (more on request); for driver operation a PC with the following minimum requirements is needed: min. Windows 7, min. 8 GB RAM, min. AMD Athlon 240GE or Intel Core i5 6200U or comparable
- Power demand < 30 W, PSU included
- BNC front connectors, 50 Ω coaxial type
- USB2.0 Type B communication connector
- DC power supply, mains connection type F (others on request)
- Factory calibration

# Scope of delivery

- Tensormeter RTM1 19" rack-mountable device
- Software package with Windows drivers and client code examples for LabVIEW and Python
- Tensormeter RTM1 User Guide incl. TCP Commands, Operation Parameter Guidelines
- USB cable, power cable, power supply 24V / 40W

Tensor Instruments is a division of HZDR Innovation GmbH, Bautzner Landstr. 400, 01328 Dresden, Germany. +49 (0) 351 260 2590 | info@hzdr-innovation.de | www.tensorinstruments.com

Tensormeter is a patented development of Helmholtz-Zentrum Dresden-Rossendorf e.V. | All rights reserved. 2021.